

High-sensitivity, Low-noise Mini Optic Fiber Spectrometer

ATP3000/ATP3040

Feature:

- Resolution: 0.01-2 nm
- Detector pixels: 2048
- High sensitivity: 1300V (lx*s)
- Response range: from UV to NIR
- Maximum readout speed: 10MHz max
- Optical path: crossed C-T
- Integration time: 1ms-65535ms
- Power supply: USB port
- Light input connector: SMA905 or free space
- Data output connector: USB2.0(High speed) or UART

Application:

- Small volume, fast spectrophotometer
- Spectral analysis/radiation spectrophotometer analysis
- Transmission, absorbance measurement
- Reflectance detection
- Ellipsometer
- Laser wavelength measurement
- UV, VIS and short wave NIR wavelength measurement

Description:

Optosky ATP3000 Series are low-noise, high-resolution optic fiber spectrometer. It employs 2048pixels CCD sensor spectral response covering UV to NIR wavelength. Readout speed reach up to 10MHz, low noise signal process circuit, and fast and accurate spectral data collection.

ATP3040 further improve SNR and dynamic range based on ATP3000, so measuring results have high stability and reliability, and they do not change with operating temperatures.

ATP3000 supplies power directly via USB port, and outputs data via USB2.0/UART, and it's convenient to apply integration.

PN	CCD pixels	Cooled or uncooled CCD
ATP3000	2048pixels	Uncooled CCD
ATP3040	4096pixels	Uncooled CCD




Product data information is current as of publication data. Products conform to specifications per the terms of Optosky Standard warranty.

1

Copyright © Optosky Photonics Inc., 2015
 Floor 15th, Bld. A04, 3rd Software Park, Jimei,
 Xiamen, 361005, China
 Tel: +86-592-6102588


Performance parameters:



Detector:		
Type		Linear CCD sensor
Spectral range		180-1100 nm
Resolution		0.01-2 nm
Effective pixels		2048pixels（ATP3000） or 4096 pixels（ATP3040P）
Pixel size		14μm×200μm （ATP3000） 或 7μm×200μm（ATP3040P）
Sensitivity		1300（lx·s）
conversion efficiency		25uV/e-
Dark noise	ATP3000	8.3 RMS @ 25℃
Optical parameters		
Wavelength ranges		180-1100 nm （available in custom wavelength）
Optical resolution		0.01-1 nm （decide on slit and spectral range）
SNR	ATP3000	>1000： 1
Dynamic range	ATP3000	10000： 1
Optical path parameters		
Optical path		f/4 crossed C-T
Confocal distance		77.5 mm for incidence / 112 mm for output
entrance slit width		5、10、25、50、100、150、200 μm (optional), available in custom width)
incident connector		SMA905 connector or free space
Electrical parameters		
Integration time		1 – 65535 ms
Output data port		USB 2.0
ADC bits depth		16 bit
Power supply		DC 5V±10%
Operating current		<350 mA（uncooled）、 <3.0A（cooled）
Storage temperature		-20℃ to +70℃
Operating temperature		-10-45℃
Operating humidity		< 90%RH
Physical parameters		
size		170×110×52 mm ³
weight		0.8 kg
Sealing		Anti-sweat

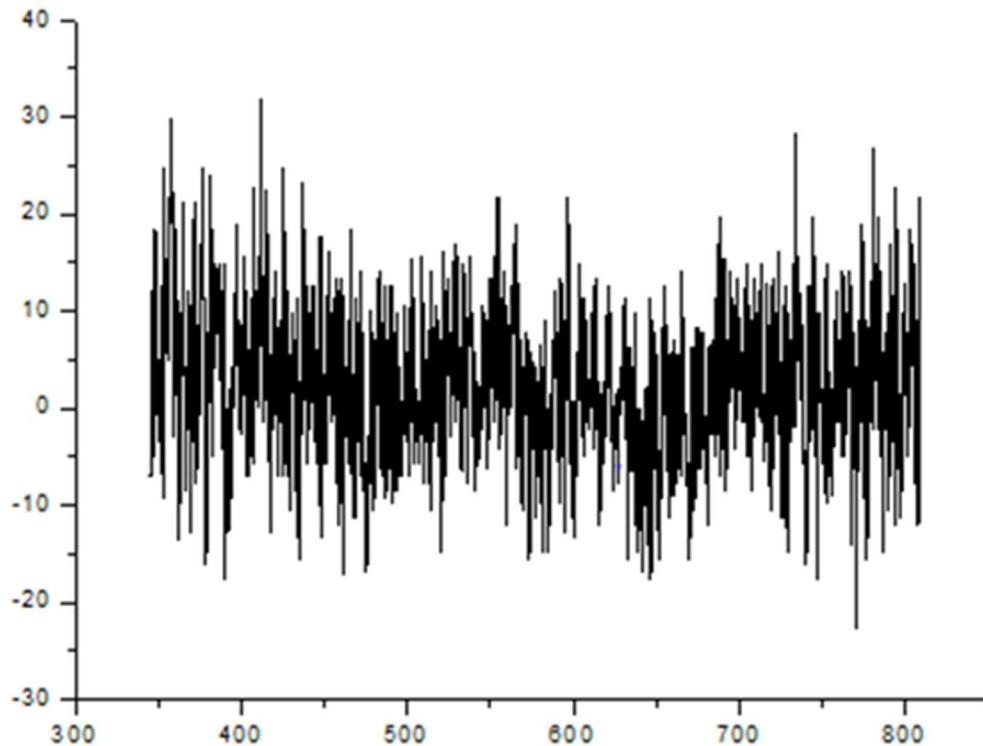

ATP3000 dark noise

Mechanical dimension:

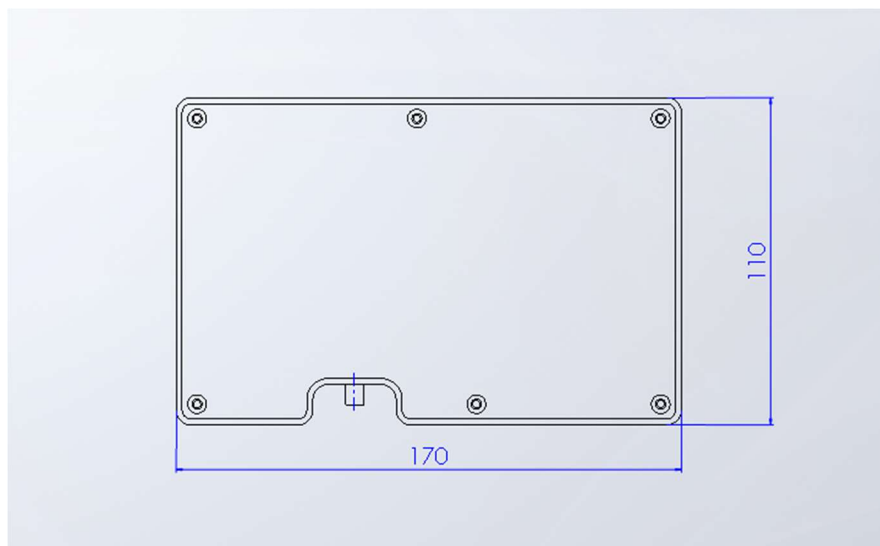

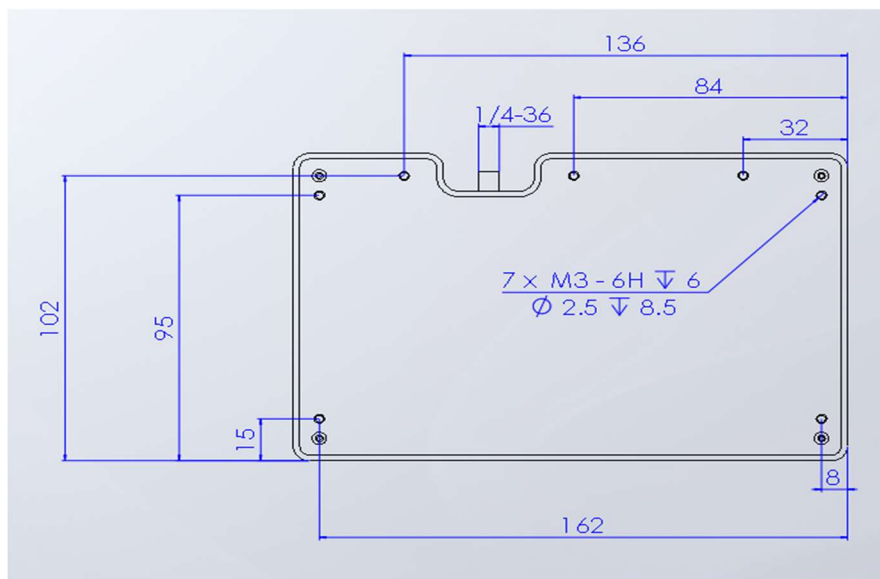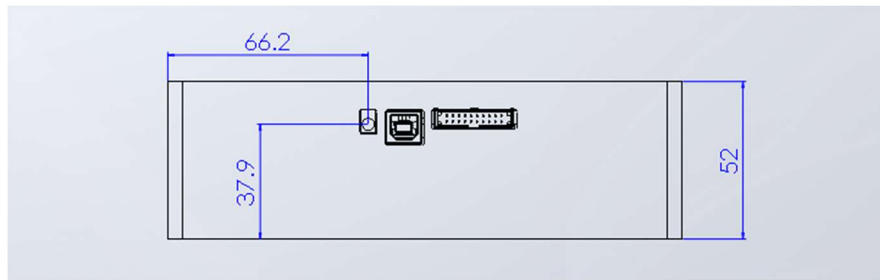

Electrical port:

Parameter	Min	Typ	Max	Unit
Power Supply				
Operating voltage range	4.5	5	5.5	V
Operating current	200	500	2000	mA
Logic Inputs(3.3V LVTTTL, Five-volt tolerant)				
High level input voltage	1.7		3.6	V
Low level input voltage	-0.3		1.0	V
Logic Output(3.3V LVTTTL)				
High level output voltage	2.4			V
Low level output voltage			0.4	V

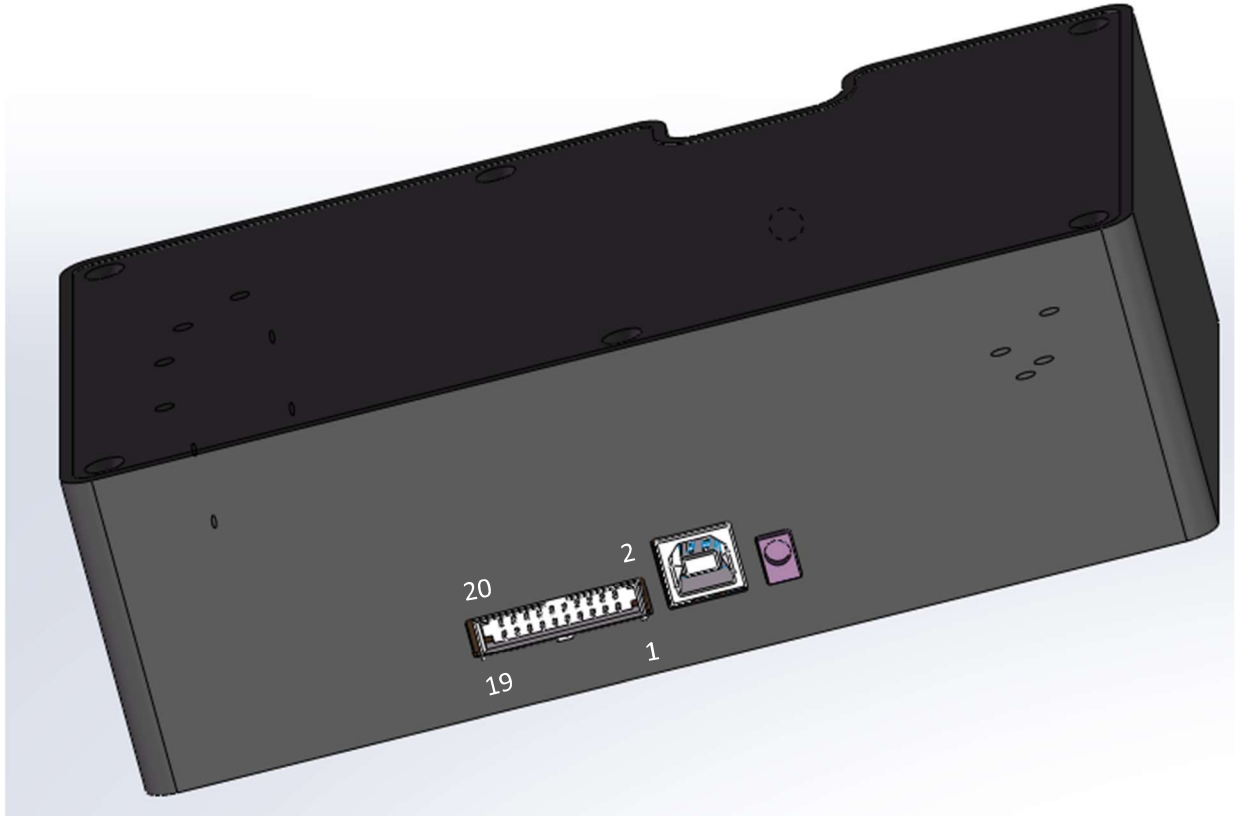


Table 1 Electrical Pin-Out

Pin#	Description	I/O	Function Description
1	VCC	/	Power Supply, $5V \pm 0.5$,
2	VCC	/	Power Supply, $5V \pm 0.5$,
3	GND	/	Ground
4	GND	/	Ground
5	LD_TX	Output	UART Transmit signal LVTTTL Logic for LD
6	LD_RX	Input	UART Receive signal LVTTTL Logic for LD
7	LD_trigger	Input	LVTTTL output trigger signal for LD
8	LD_EN	Output	LVTTTL output enable signal.for LD
9	NC	/	
10	NC	/	
11	GPIO0	Input /Output	General Purpose Software Programmable Digital Inputs/Outputs, LVTTTL Logic.
12	GPIO1	Input	General Purpose Software Programmable Digital

		/Output	Inputs/Outputs, LVTTTL Logic.
13	GPIO2	Input /Output	General Purpose Software Programmable Digital Inputs/Outputs, LVTTTL Logic.
14	GPIO3	Input /Output	General Purpose Software Programmable Digital Inputs/Outputs, LVTTTL Logic.
15	GPIO4	Input /Output	General Purpose Software Programmable Digital Inputs/Outputs, LVTTTL Logic.
16	GPIO5	Input /Output	General Purpose Software Programmable Digital Inputs/Outputs, LVTTTL Logic.
17	VCC	/	3.3V Power Output
18	GND	/	Ground
19	EXT_TX	Output	EXT UART Transmit signal LVTTTL Logic
20	EXT_RX	Input	EXT UART Receive signal LVTTTL Logic